



## Exporting Results from Tables in Empower

Create a new Export method: Browse projects->File->New Method->Export Method. Fill the information in the FIELDS tab at the right places: Check the Table->Export Table Data and uncheck the “Export Row Numbers”. Edit the Table as needed. Add “Sample Information” as needed.

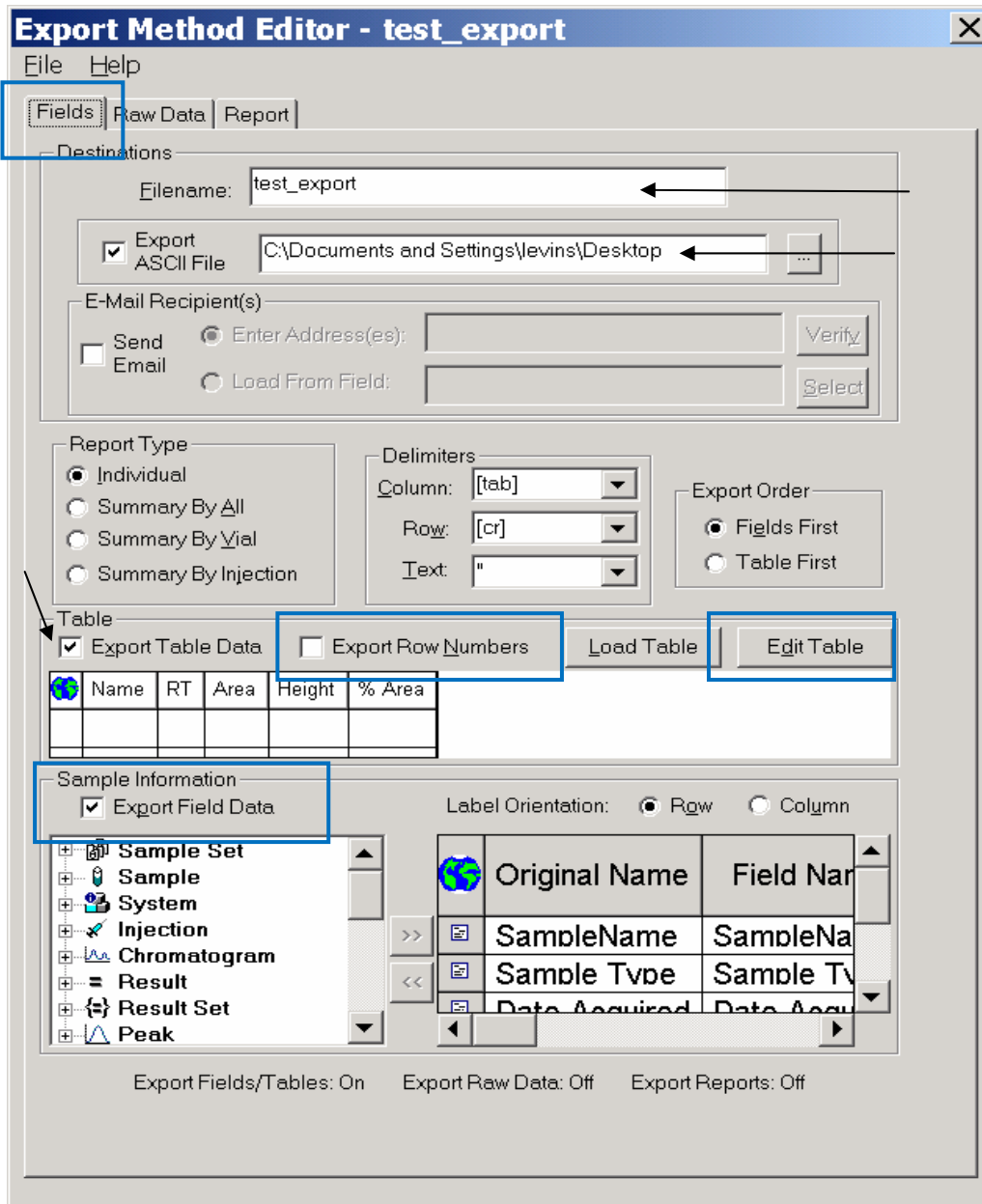


Save the method and go to your results. Select the results to be exported and Right-Click->Export. The following screen will be opened: select the Export Method that you just created.


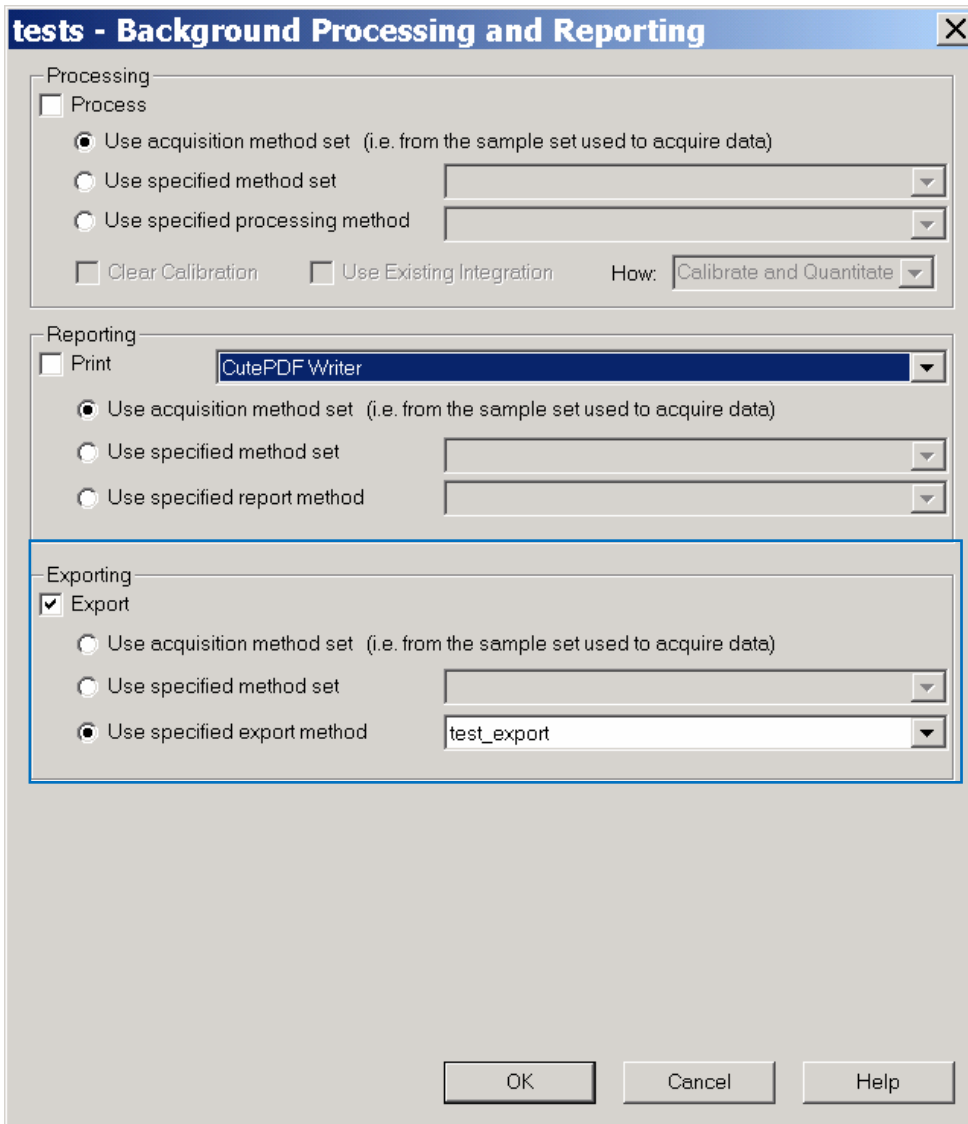


A tabulated text file by the name you entered in the *Export Method* will appear at the address selected in the *Export Method: name.txt*


Open Excel->File Open (Select: All Files).

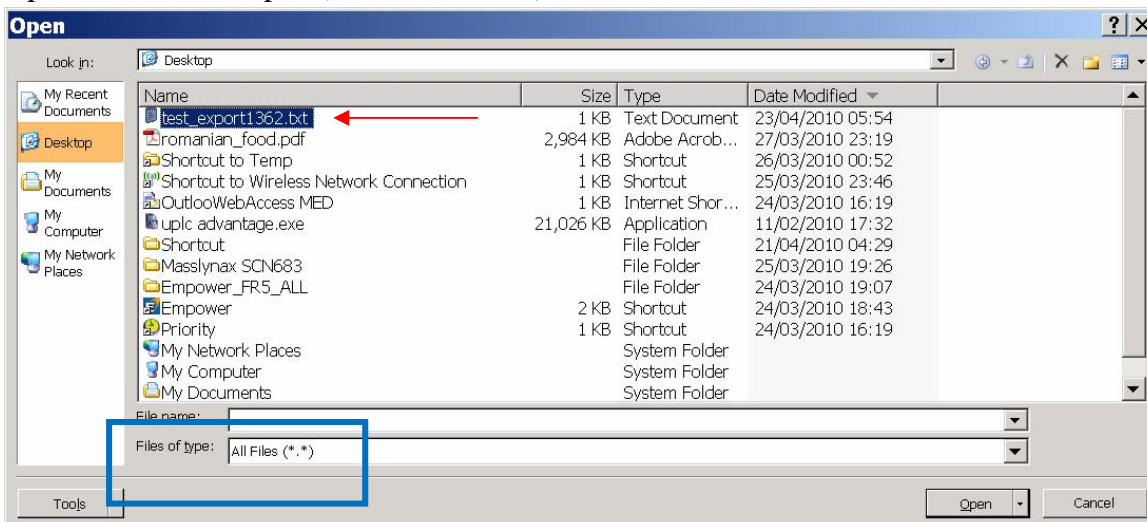

A wizard will be opened in which you do not need to change anything, just go NEXT in the following 3 steps:

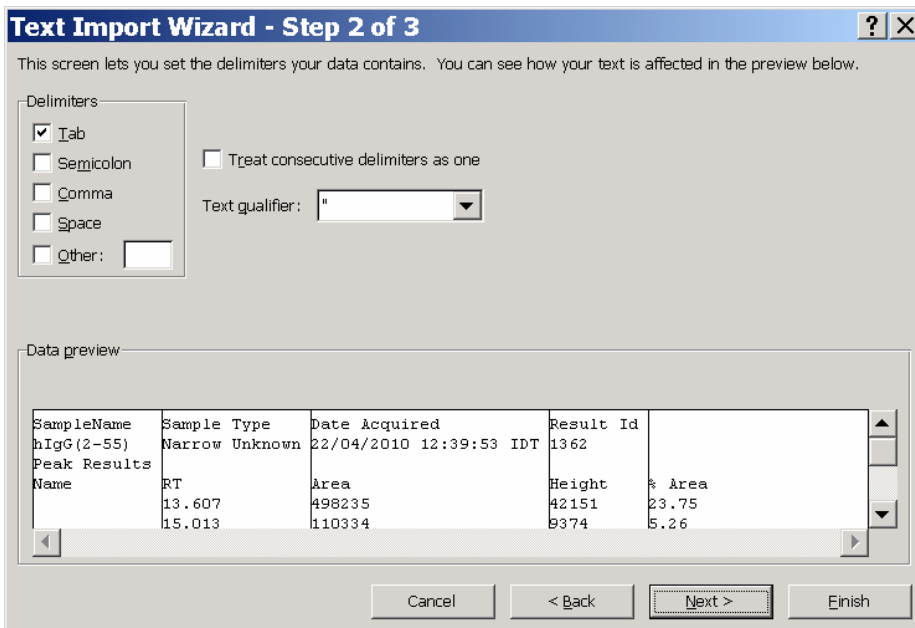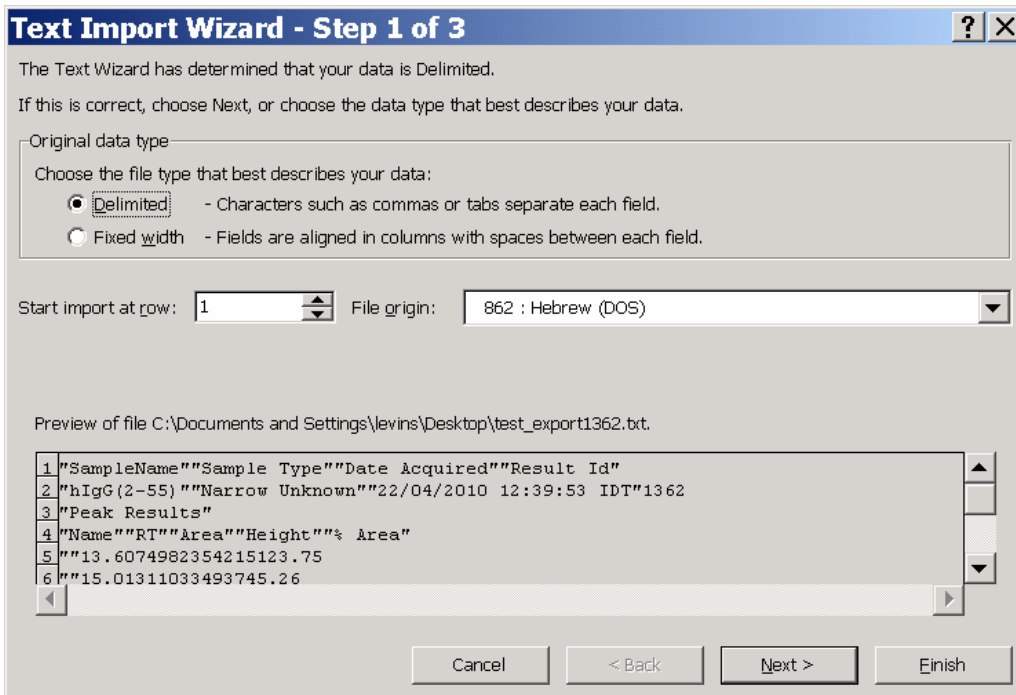

Click on Finish in the last screen:

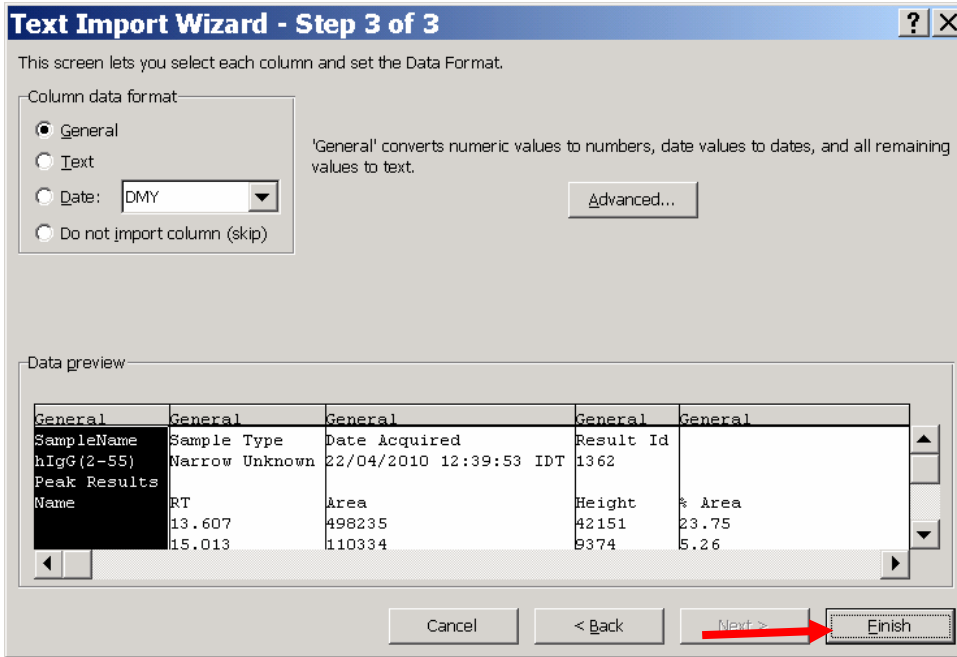

The data will appear in Excel ready and tabulated properly

	A1		fx	SampleName		
	A	B	C	D	E	F
1	SampleName	Sample Type	Date Acqu	Result Id		
2	hIgG(2-55)	Narrow Ur	22/04/201	1362		
3	Peak Results					
4	Name	RT	Area	Height	% Area	
5		13.607	498235	42151	23.75	
6		15.013	110334	9374	5.26	
7		16.483	515840	41509	24.59	
8		16.901	257752	19092	12.29	
9		17.567	102210	7505	4.87	
10		18.221	18063	1539	0.86	
11		19.617	392963	24497	18.74	
12		20.374	33628	1827	1.6	
13		21.829	37772	2029	1.8	
14		24.189	130603	5247	6.23	
15						
16						